Long-arm Biotin Conjugation Kit

Product Information

Cat No	Product Name	Size
BIOS-2022K	Long-arm Biotin Conjugation Kit	1 mg

Description

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Note

1. Store the kit at 2-8℃, do not freeze.
2. The components of the kit may be upside down during transportation, causing liquid or dry powder reagents to stick to the tube wall or bottle cap. Please centrifuge before use to make the liquid or dry powder reagent attached to the tube wall or bottle cap settle to the bottom of the tube.
3. Biotin needs to be prepared for immediate use, and the dissolved Biotin cannot be stored for a long time.

4. The buffer in the purification column contains the toxic component sodium azide (NaN_3). Avoid contact with skin, eyes and mucous membranes when using it.
5. DMSO is slightly toxic, permeable to human skin and irritating to the eyes. Avoid contact with skin, eyes and mucous membranes when using it.
6. The dialysis, concentration and concentration determination of the antibody before labeling will all cause the loss of the amount of antibody. Therefore, when preparing the antibody before labeling, consider the most appropriate amount of antibody according to the specific situation.

Contact Us

QQ:499854788

3494243873

WeChat: 13681256816

17731100244

Email: info@biotyscience.com

Tel: 400-669-8850